



Paroxetine and Desfluoro Analogue

ACE[®]
Ultra-inert
UHPLC & HPLC Columns


Application #AN3890

Conditions


Column: ACE 5 CN
Dimensions: 150 x 4.6 mm
Part Number: ACE-124-1546
Mobile Phase: 20 mM ammonium formate pH 3.0/MeOH (60:40 v/v)
Flow Rate: 2 mL/min
Injection: 20 µL
Temperature: Ambient
Detection: UV, 295 nm

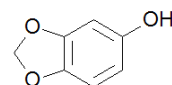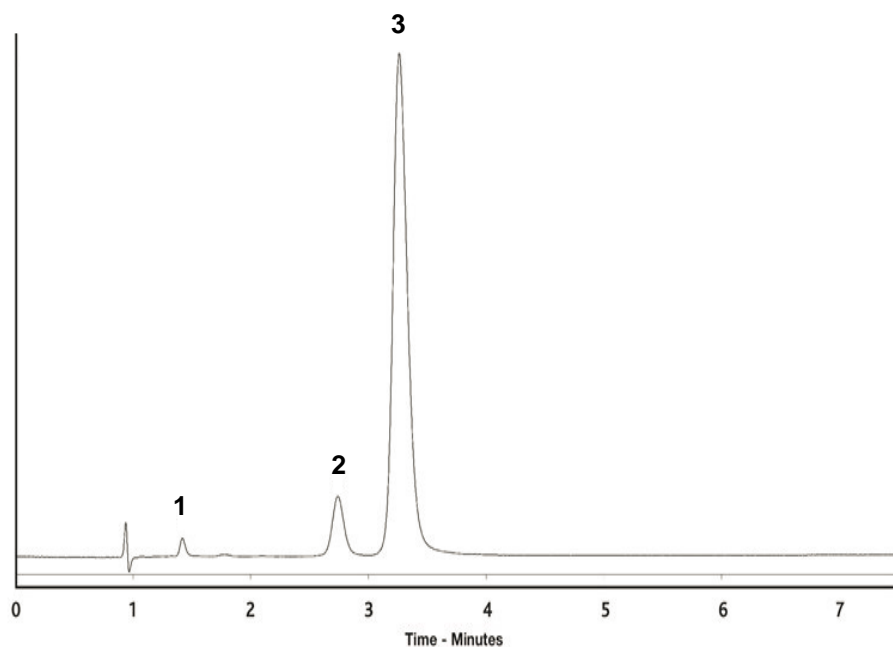

1. Sesamol

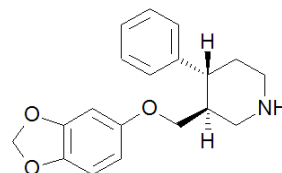

2. Desfluoroparoxetine

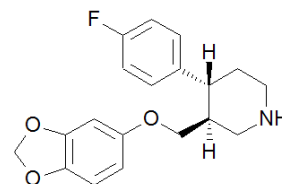

3. Paroxetine

Reproduced with permission of School of Pharmacy, University of Sunderland, UK
© 2015 ACT Ltd. All rights reserved. All products are available worldwide.

For more information contact your local ACE distributor or visit
www.ace-hplc.com or email: info@ace-hplc.com


ACE[®]
UHPLC & HPLC
Columns