



Conditions


| | |
|---------------|--|
| Column: | ACE Excel 5 NH2 |
| Dimensions: | 250 x 4.6 mm |
| Part Number: | EXL-1214-2546U |
| Mobile Phase: | MeCN:H ₂ O 80:20 v/v |
| Flow Rate: | 1.5 mL/min |
| Temperature: | 35 °C |
| Detection: | Refractive Index |
| Sample: | Carbo Energy Drink and Sugar standards |

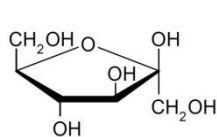


1. Fructose

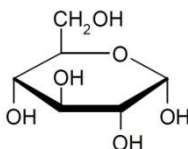


2. Glucose

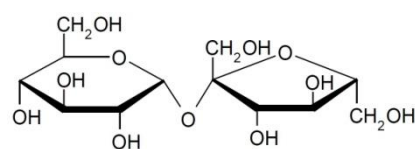


3. Sucrose

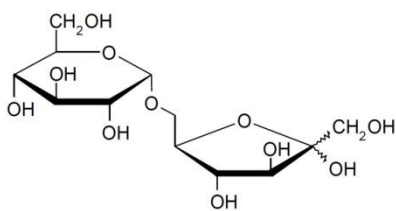

4. Palatinose

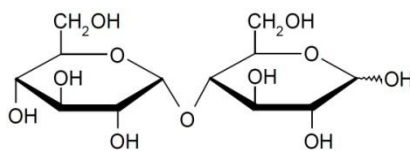


5. Maltose

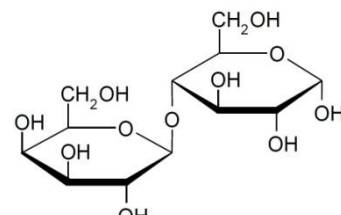


6. Lactose

Blue Exotic Carbo Energy Drink

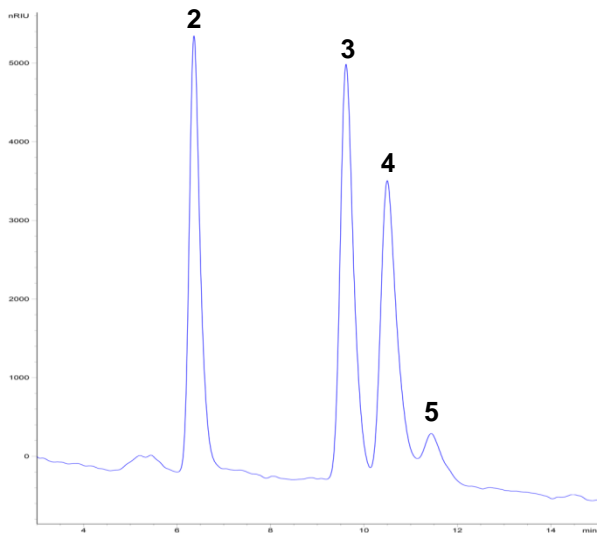

Sugar standards

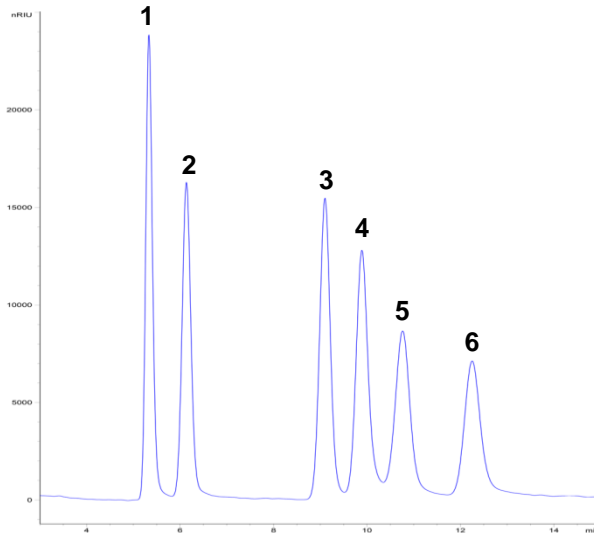

Reproduced with permission of Clinical Nutrition, S.A.U., Barcelona, Spain

© 2016 ACT Ltd. All rights reserved. All products are available worldwide.